



73 Mall Drive • Commack, NY 11725 • www.accu-scope.com • Phone (631) 864-1000 • Fax (631) 543-8900

3032 Metal Halide Collimating Adapter

ASSEMBLY INSTRUCTIONS

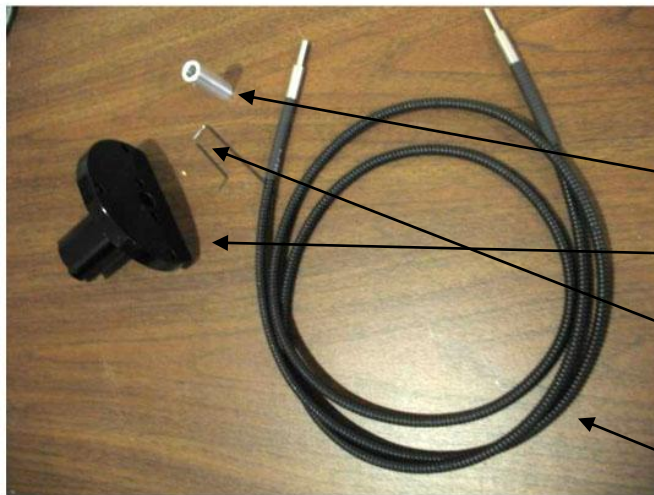


FIG. 1

Components

(FIG. 1)

Silver ferrule

Collimating adapter
(half moon plate & attached lens tube)

.005 Allen wrench

1/16 inch Allen Wrench (*not shown*)

Light guide



FIG. 2

Assembly

(FIG. 2)

Take the silver ferrule and place it on the light guide tip.



FIG. 3

Assembly *(continued)*

(FIG. 3 & 4)

Align the tip of the light guide so that it is flush with the end surface of the ferrule.

On the trailing edge of the ferrule there is a small set screw that tightens with the .005" Allen wrench.

Tighten the ferrule to the light guide.

Locate the other set screw in a hole on the left side of the lens tube. It is approximately 1 inch from the surface of the half moon plate (FIG. 4).



FIG. 4

(FIG. 5)

NOTE: BEFORE you install the collimating adapter, insert the silver ferrule into the lens tube until it stops. The trailing edge of the ferrule should end up about 1/4" inside the half moon plate.

Using a 1/16" Allen wrench, tighten the set screw that is on the left side of the lens tube. This set screw secures the light guide at the proper position.

(FIG. 6)

Assembly complete. Attach the dovetail slider and install the collimating adapter to the 3032.



FIG. 5



FIG. 6