



## **3032 Stage Replacement**

1. Disconnect the lamp power cord from the rear of the microscope.
2. Remove the four screws securing the lamp post to the stage. **NOTE:** There may be one or more shims between the lamp post and the stage. If so, reuse them in the same location(s) when installing the lamp post onto the new stage.
3. Remove the screw-head covers over the four screws securing the stage to the microscope frame. Remove the screws and lift the stage off the frame.
4. Install the lamp post onto the stage.
5. Install the stage onto the microscope frame and install the new screw-head covers.
6. Reconnect the lamp power cord to the microscope base.