



# EXC-500 Microscope Series

## 2x Diffuser Supplement

*2x Diffuser (Swing-in, bayonet mount) for the EXC500 Microscope, Part # 500-3222-2x  
This component may be required with some combinations of video cameras, c-mounts and monitors.*



This component is comprised of two flanges; a semi-permanent mounting flange and a swing-in diffuser. The lower flange has a four-tooth bayonet arrangement. Placing it on the light port as shown below will allow it to seat down onto the light port.



Next the assembly is rotated 45 degrees. This will engage the bayonet teeth. When correctly installed, the thumb screw on the lower flange should face 3 o'clock. Lightly tighten the thumb screw.



To temporarily remove the diffuser from the light path swing the upper flange CCW.