

EXC-500 Microscope Series

Objective Installation Sequence



To install the objectives:

1. Turn the coarse adjustment knob until the revolving nosepiece is at its lowest position.
2. Remove the nosepiece cap closest to the position of the RED DOT.
3. Beginning with the lowest magnification objective thread the objective onto the nosepiece opening, then rotate the nosepiece clockwise and thread the other objectives from low to high magnification.
4. You can remove the red dot once all the objectives are installed.

NOTE

- Always rotate the nosepiece by using the knurled nosepiece ring.
- Keep the covers on any unused nosepiece openings to prevent dust and dirt from getting inside.